


700 Series Triple Beam Balances, Fig 1


Click on the key number. The part information will be on the next page shown

Click here for items not shown on the drawing

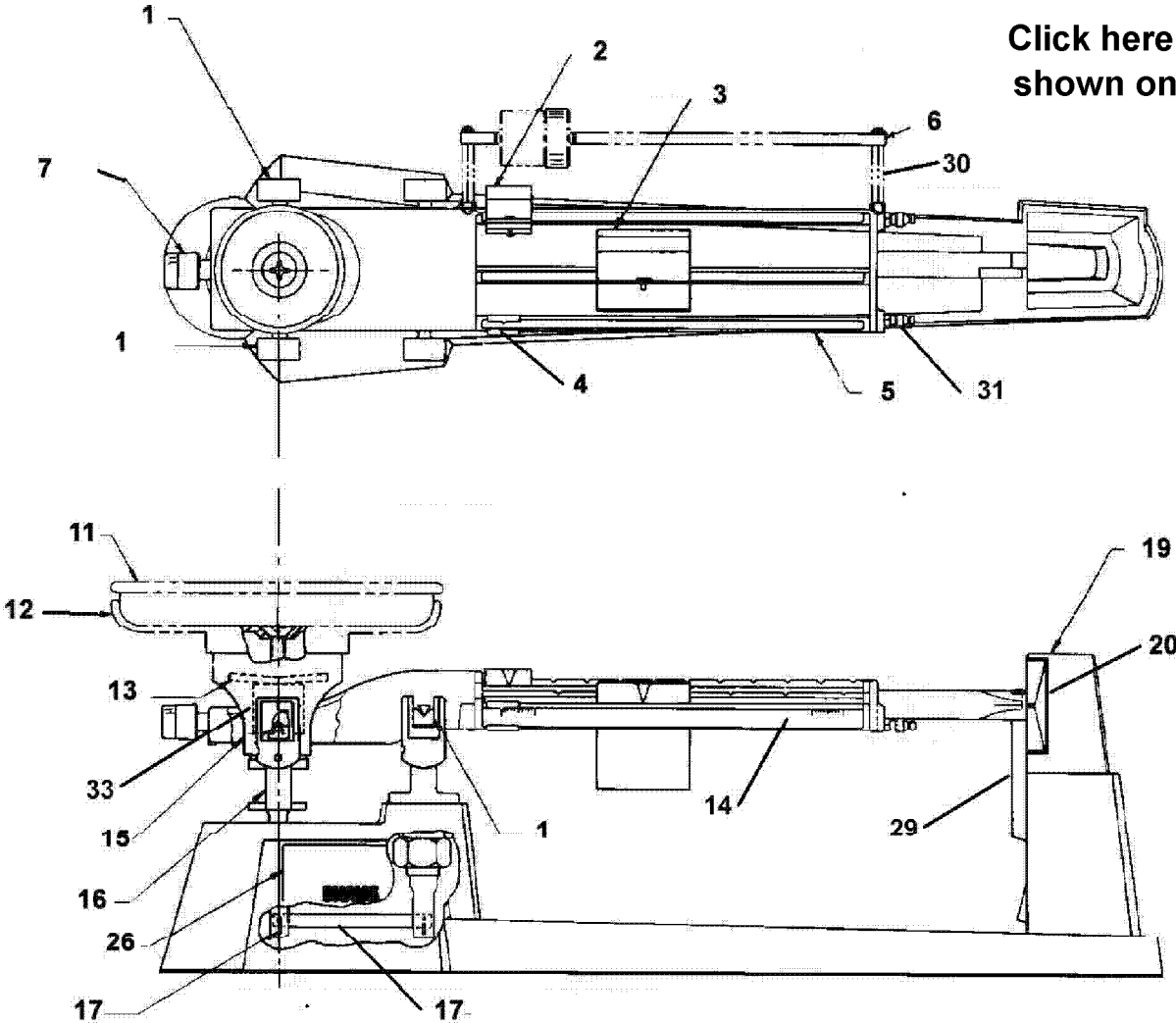


700 Series Triple Beam Balances, Fig. 2


Click on the key number. The part information will be on the next page shown

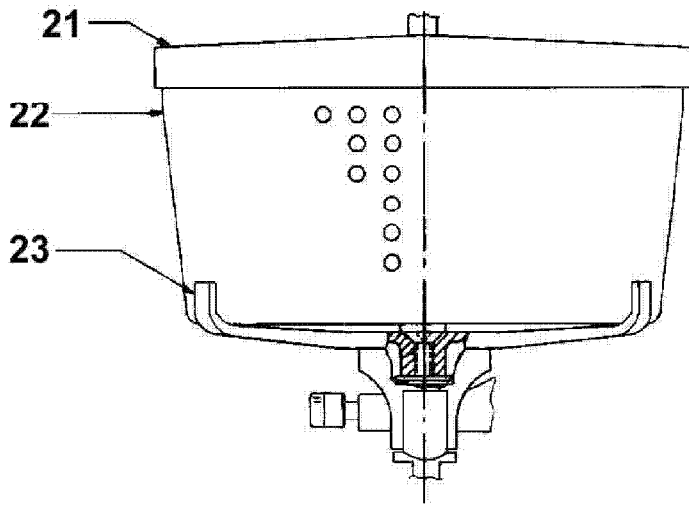


Model 730-00

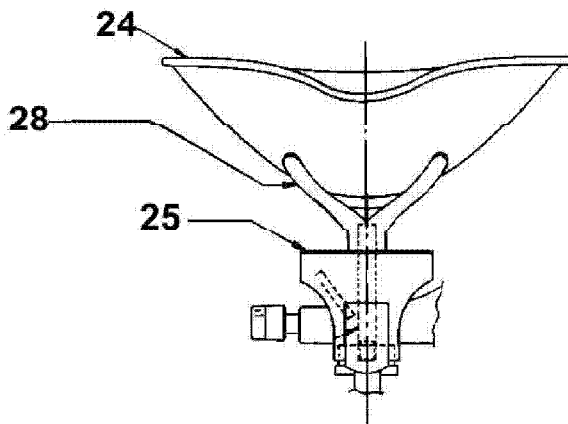

Model 720-00, 720-S0

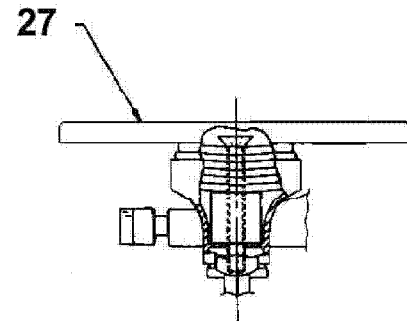


Model 760-00

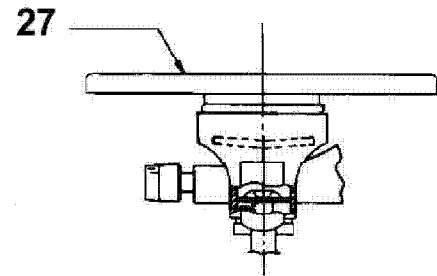

Model 750-00

1600 Series Dial balances, Fig 3.

Click on the key number. The part information will be on the next page shown

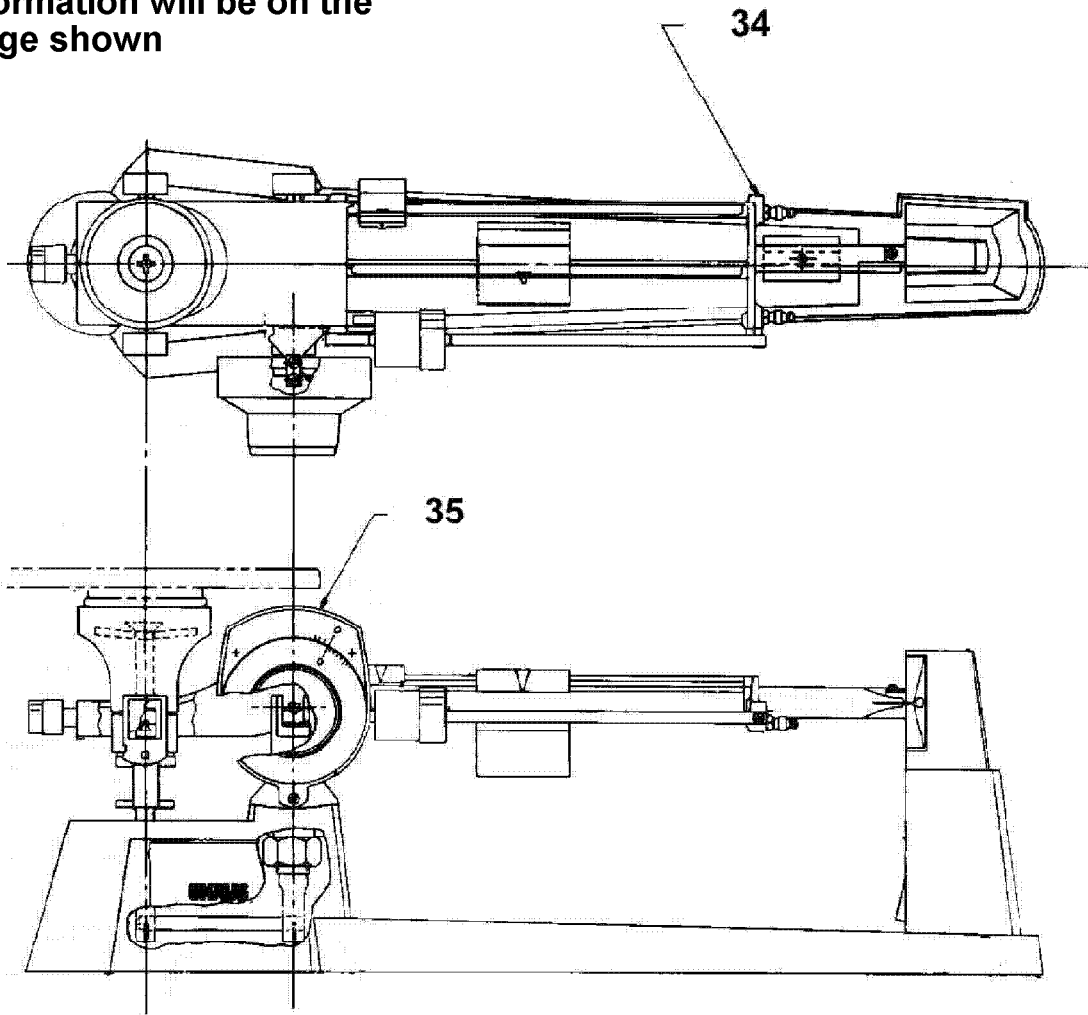



AfterMarket
Parts • Service • Support • Accessories

ServiceParts List

Series: Triple Beam Balance

Status: Active

Model: 700, 800

Key	Figure	Part No.	Description
01	01	80251252	Bearing and Cover Kit, TBB
02	01	80251328	Poise Assy, 100g, TBB, DOG 1977 on
03	01	80251327	Poise Assy, 500g, TBB, DOG
04	01	80251326	Poise Assy, 10g, TBB
05	01	3266-30	Beam Assembly, 700 series, 10g poise with Tare
06	01	80251277	Tare Beam Assy, TBB, DOG
07	01	80251250	Balance Compensator Kit, TBB, HVD
11	01	80251313	SS Pan, TBB, HVD
12	01	80251317	Cross, TBB, HVD
13	01	1097-03	Balancing Slug, No Hole 700, 1400,310, 311
14	01	80251255	Graduated Inserts Kit, TBB
15	01	80251310	Balance Cup, TBB, HVD
16	01	80251311	Parallel Loop Assembly, TBB, HVD
17	01	80251262	Check Rod Kit, TBB
19	01	80251264	Trig Loop Assy with Dial Plate and Magnets, TBB
20	01	80251266	Dial Plate, TBB
26	01	80251278	Label, TBB
26	01	80251279	Label, TBB, DOG
29	01	80251258	Damper Vane Kit, TBB
30	01	80251275	Tare Beam lug TBB, HVD
31	01	80251261	Suspension Pivots Kit, TBB



AfterMarket
Parts • Service • Support • Accessories

ServiceParts List

Series: Triple Beam Balance

Status: Active

Model: 700, 800

Key	Figure	Part No.	Description
21	02	3052-00	Specimen Box Cover
22	02	3043-20	Specimen Box without cover
23	02	2570-21	Cross Assembly
24	02	80250400	(Sold as Accessory) Stainless Steel Scoop without Foot 700 1400
24	02	1101-20	(Sold as Accessory) Scoop, Polypropylene
25	02	80251318	Balance Cup Cover and Stud, TBB
27	02	80251314	Scale Plate Assembly, TB, HVD
28	02	80251333	Fork, TBB



AfterMarket
Parts • Service • Support • Accessories

ServiceParts List

Series: Triple Beam Balance

Status: Active

Model: 700, 800

Key	Figure	Part No.	Description
34	03	3266-40	Beam Assembly without Poises, 1600
35	03	80251323	Dial Assembly, Complete (DOG only)



AfterMarket
Parts • Service • Support • Accessories

ServiceParts List

Series: Triple Beam Balance

Model: 700, 800

Status: Active

Key	Figure	Part No.	Description
	NS	701-00	500g Attachment Weight, 700
	NS	80251293	Shipping Kit, Complete, DOG
	NS	80252228	Shipping Kit, Complete, TBB
	NS	3258-00	Manual, Instru. Triple, EN
	NS	80251271	Packing washer, TBB, HVD
	NS	9751-18	Bag of Shot
	NS	702-00	1000g Attachment Weight, 700